

# Extraction of Aflatoxins and Ochratoxin from Dried Chili Using ISOLUTE® Myco Prior to LC-MS/MS Analysis

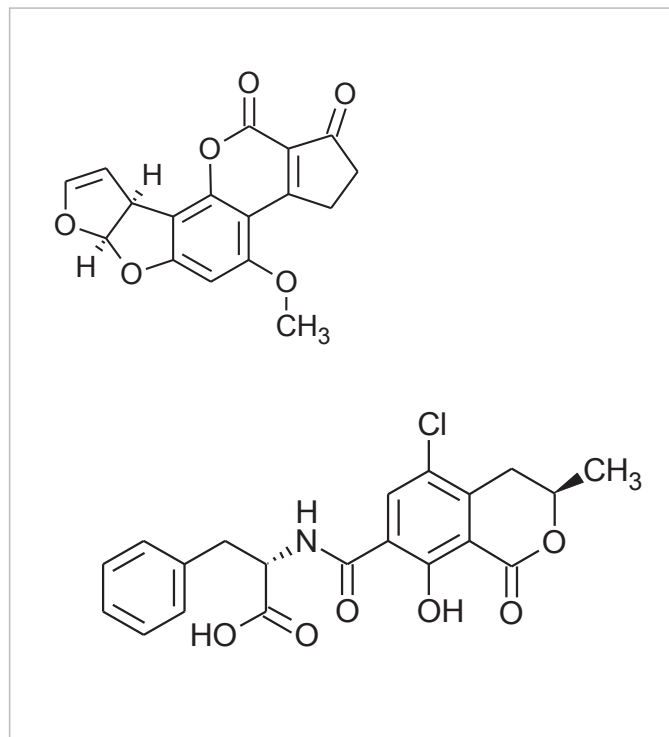


Figure 1. Structures of Aflatoxin B1 and Ochratoxin A

## Introduction

This application note describes a solid phase extraction (SPE) protocol for the extraction of a range of mycotoxins from dried chili (pimiento) using ISOLUTE® Myco with LC-MS/MS analysis.

Mycotoxins are toxic metabolites produced by fungal molds on food crops. Regulation and legislation for testing of mycotoxin contamination has established which mycotoxins are prevalent on a wide variety of food crops. This application note describes an SPE protocol appropriate for LC-MS/MS analysis of a range of mycotoxins found on chili.

The method described in this application note achieves high recoveries of relevant mycotoxins from dried chili (pimiento) with %RSDs and LOQs that meet the requirements set in European Union regulations for measurement of these analytes. ISOLUTE® Myco solid phase extraction columns provide robust, reliable sample preparation for multiple mycotoxin classes from a wide range of foodstuffs. Using a single, easy to use sample preparation product, along with optimized matrix specific application notes, scientists can prepare diverse food/crop samples for analysis by LC-MS/MS.

## Analytes

Aflatoxin B1, aflatoxin B2, aflatoxin G1, aflatoxin G2, ochratoxin A

## Sample Preparation Procedure

**Format:** ISOLUTE® Myco 60 mg/3 mL Columns (Tabless), part number 150-0006-BG

### Sample Pre-treatment

- Sample processing:** Grind the sample (50 g) with a burr-grinder or equivalent device. Store ground sample in a sealed container at room temperature until required.
- Extraction:** Mix the ground sample (5g) with 80% acetonitrile (aq) (20 mL). Place the sample pre-treatment tube on a shaking table for 30 mins. Transfer the extract to a 50 mL centrifuge tube and centrifuge at 4000 g for 10 minutes.
- Dilution:** Take the supernatant (2 mL), transfer to a new 50 mL centrifuge tube and dilute with water (32 mL). Centrifuge diluted extract at 4000 g for a further 10 minutes.

## Solid Phase Extraction

Use flow rates of 1 mL min<sup>-1</sup> throughout

<b>Condition:</b>	Condition the column with acetonitrile (2 mL)
<b>Equilibration:</b>	Equilibrate column with water (2 mL)
<b>Sample Loading:</b>	Load pre-treated sample (3 mL) onto the column at a maximum flow rate of 1 mL min <sup>-1</sup> (gravity load is recommended)
<b>Interference Wash 1:</b>	Wash the column with water (2 x 2.5 mL)
<b>Interference Wash 2:</b>	Wash the column with 10% acetonitrile (aq) (2 x 2.5 mL)
<b>Drying:</b>	Dry the column for 30 seconds at maximum vacuum, -0.5 bar/7 psi
<b>Elution 1:</b>	Elute with 0.1% formic acid in 40% acetonitrile (aq) (2 mL)
<b>Elution 2:</b>	Elute with 1.0% ammonia (conc.) in methanol (2 mL)
<b>Post Elution:</b>	Dry the eluate in a stream of air or nitrogen using a SPE Dry (35 °C, 20 to 40 L min <sup>-1</sup> ) or TurboVap® LV (15 bar at 35 °C for 40 min). Reconstitute in 0.1 % acetic acid in acetonitrile : methanol : water (1 : 1 : 8, v/v/v, 1mL). Syringe-filter using a 0.2 µm PTFE membrane prior to analysis.

## HPLC Conditions

<b>Instrument:</b>	Shimadzu Nexera UHPLC (Shimadzu Europe GmbH)
<b>Column:</b>	Kinetex XB-C18 50 x 2.1 mm 2.6 µm dp (Phenomenex, Macclesfield UK)
<b>Mobile Phase:</b>	A: 1 mM ammonium acetate, 0.5% acetic acid B: 1mM ammonium acetate, 0.5% acetic acid in 95% methanol (aq)
<b>Flow Rate:</b>	0.45 mL min <sup>-1</sup>
<b>Injection:</b>	20 µL
<b>Gradient:</b>	Initial 20 % B, hold 1.0 min linear ramp to 73 % B in 6 min linear ramp to 100 % B in 0.2 min, hold 2.3 min linear ramp to initial conditions in 0.2 min hold 2.3 min, total run time 10.0 min
<b>Column temperature:</b>	40 °C
<b>Sample temperature:</b>	15 °C

**Table 1.** Typical retention times for a range of mycotoxins using the LC-MS/MS method described.

Compound	Retention time (min)
aflatoxin G2	3.3
aflatoxin G1	3.6
aflatoxin B2	3.9
aflatoxin B1	4.1
ochratoxin A	6.1


## MS Conditions

Ions were selected in order to achieve maximum sensitivity, and the MS was operated in positive ion polarity mode, using multiple reaction monitoring.

<b>Instrument:</b>	AB Sciex Triple Quad 5500 (Warrington, UK)
<b>Source:</b>	Turbo-V ESI
<b>Desolvation temp.:</b>	500 °C
<b>Curtain gas:</b>	30 psi
<b>Spray voltage:</b>	+5.0 kV
<b>Gas 1:</b>	60 psi
<b>Gas 2:</b>	60 psi
<b>Collision gas:</b>	7 psi

**Table 2.** Positive Ion Mode - MRM Parameters

MRM transition	RT	Compound ID	DP, V	EP, V	CE, V	CXP, V
331.1>313.1	3.3	aflatoxin G2 1	100	10	33	12
331.1>245.1	3.3	aflatoxin G2 2	100	10	41	12
331.1>257.1	3.3	aflatoxin G2 3	100	10	41	12
329>243.1	3.6	aflatoxin G1 1	80	10	37	12
329>200	3.6	aflatoxin G1 2	80	10	53	12
315.1>287	3.6	aflatoxin B2 1	100	10	35	12
315.1>259.1	3.9	aflatoxin B2 2	100	10	40	12
315.1>243.1	3.9	aflatoxin B2 3	100	10	51	12
313.1>285	4.1	aflatoxin B1 1	100	10	31	18
313.1>241.1	4.1	aflatoxin B1 2	100	10	49	18
313.1>185	4.1	aflatoxin B1 3	100	10	65	18
404.1>239	6.1	ochratoxin A 1	165	10	32	12
404.1>221	6.1	ochratoxin A 2	165	10	47	12
404.1>102	6.1	ochratoxin A 3	165	10	84	12

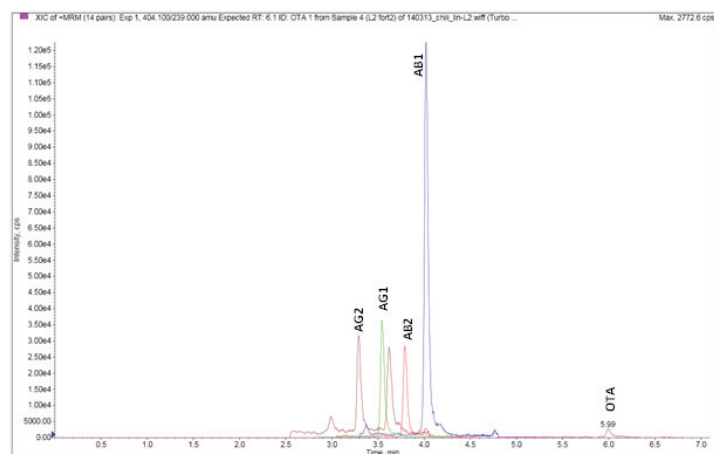

**Figure 2.** Extracted ion chromatogram (positive ion mode) using ISOLUTE® Myco protocol at 5 µg kg<sup>-1</sup> (aflatoxins and ochratoxin A) from chili

## Validation Criteria

Method linearity was determined using matrix-matched calibration standards in six replicates over six levels; the ranges are shown below.

Analytes	Working Range, $\mu\text{g kg}^{-1}$ ( $\mu\text{g } \mu\text{L}^{-1}$ on column)
aflatoxin B1, aflatoxin B2, aflatoxin G1, aflatoxin G2, ochratoxin A	0.5 to 80 (0.046875 to 7.5)
	5.0 to 80 (0.46875 to 7.5)


LOQ was determined from the lowest matrix-matched standard meeting EU repeatability and recovery criteria; or estimated from the S/N ratio of the blank where incurred levels of mycotoxins were present in the sample (e.g AB1).

Repeatability (%RSD<sub>r</sub>) was determined from single acquisitions of 5 SPE replicates of a single sample extraction. The RSDs generated gave close agreement when a single sample was extracted.


Recovery was determined as a % of ISOLUTE® Myco extract spike before sample prep to spike after at the EU MRL.

## Results

The extracted ion chromatogram in figure 2 demonstrates chromatography at  $5 \mu\text{g kg}^{-1}$  (aflatoxins and ochratoxin A) from a spiked extraction of 2 g ground chili. Good linearity was achieved for all analytes in all both matrices as demonstrated in the example charts shown in **Figures 3 and 4**.

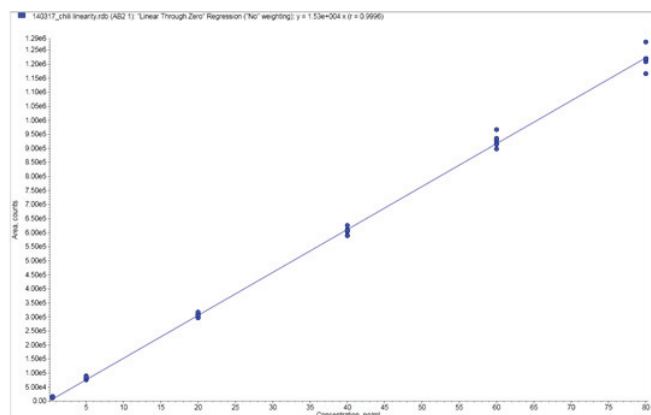

**Figure 3.** Calibration curve for aflatoxin B2 from ground chili using the ISOLUTE® Myco protocol from 0.5 – 80  $\mu\text{g kg}^{-1}$

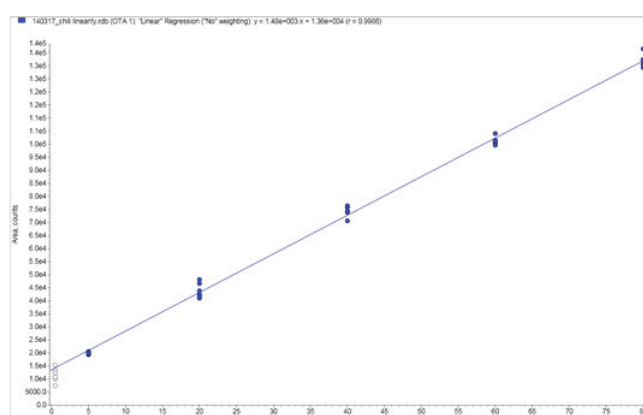

**Figure 4.** Calibration curve for ochratoxin A from ground chili using the ISOLUTE Myco protocol from 5 – 80  $\mu\text{g kg}^{-1}$

All analytes extracted using the ISOLUTE Myco protocol achieved the limits of quantities and recovery required by the current European standards for mycotoxin analysis as shown in **Table 3**.

**Table 3.** Analyte recovery and limit of quantitation data for a range of mycotoxins from chili using the ISOLUTE® Myco protocol

Analyte	$r^2$	LOQ / $\mu\text{g kg}^{-1}$		%RSD <sub>r</sub>		Recovery %	
		Target	Actual	Target	Actual	Target	Actual
aflatoxin B1	0.9988	5	2.0	20	3.2	70 to 110	108
aflatoxin B2	0.9992	5	2.0	20	3.2	70 to 110	106
aflatoxin G1	0.9986	5	2.0	20	2.9	70 to 110	106
aflatoxin G2	0.9980	5	4.0	20	2.7	70 to 110	89.1
ochratoxin A	0.9972	15	10	20	2.7	70 to 110	108

## Alternative Sample Pre-Treatment Strategy

Weigh the ground sample (2 x 1 g) into separate 7 mL Omni International sample tubes containing ceramic 2.8 mm bead lysing matrix (approximately 2 g / n=23) and add 80% acetonitrile (aq) (4 mL). Homogenize the sample in an Omni Bead Ruptor 24 using the following settings: disruption at 5.5 m s<sup>-1</sup> for 2 x 45 s with a 10 s pause.

Centrifuge the sample tubes at 4000 x g for 10 minutes. Transfer the supernatant (2 x 2 mL) to a new centrifuge tube and dilute with water (16 mL). Centrifuge the tube at 4000 x g for 10 minutes. Proceed to solid phase extraction (page 2).

## Ordering Information

Part Number	Description	Quantity
150-0006-BG	ISOLUTE® Myco 60 mg/3 mL column (Tabless)	50
121-1016	Biotage® VacMaster™-10 Sample Processing Manifold complete with 16 mm collection rack	1
121-2016	Biotage® VacMaster™-20 Sample Processing Manifold complete with 16 mm collection rack	1
C103198	TurboVap® LV, 100/120V	1
C103199	TurboVap® LV, 220/240V	1

For the latest application notes visit [www.biotage.com](http://www.biotage.com)

### EUROPE

Main Office: +46 18 565900  
Toll Free: +800 18 565710  
Fax: +46 18 591922  
Order Tel: +46 18 565710  
Order Fax: +46 18 565705  
[order@biotage.com](mailto:order@biotage.com)  
Support Tel: +46 18 56 59 11  
Support Fax: +46 18 56 57 11  
[eu-1-pointsupport@biotage.com](mailto:eu-1-pointsupport@biotage.com)

### NORTH & LATIN AMERICA

Main Office: +1 704 654 4900  
Toll Free: +1 800 446 4752  
Fax: +1 704 654 4917  
Order Tel: +1 704 654 4900  
Order Fax: +1 434 296 8217  
[ordermailbox@biotage.com](mailto:ordermailbox@biotage.com)  
Support Tel: +1 800 446 4752  
Outside US: +1 704 654 4900  
[us-1-pointsupport@biotage.com](mailto:us-1-pointsupport@biotage.com)

### JAPAN

Tel: +81 3 5627 3123  
Fax: +81 3 5627 3121  
[jp\\_order@biotage.com](mailto:jp_order@biotage.com)  
[jp-1-pointsupport@biotage.com](mailto:jp-1-pointsupport@biotage.com)

### CHINA

Tel: +86 21 2898 6655  
Fax: +86 21 2898 6153  
[cn\\_order@biotage.com](mailto:cn_order@biotage.com)  
[cn-1-pointsupport@biotage.com](mailto:cn-1-pointsupport@biotage.com)

To locate a distributor,  
please visit our website at  
[www.biotage.com](http://www.biotage.com)

### Part Number: AN785

© 2014 Biotage. All rights reserved. No material may be reproduced or published without the written permission of Biotage. Information in this document is subject to change without notice and does not represent any commitment from Biotage. E&OE. Product and company names mentioned herein may be trademarks or registered trademarks and/or service marks of their respective owners, and are used only for explanation and to the owners' benefit, without intent to infringe.

For more information visit [www.biotage.com](http://www.biotage.com).